

MasterAmp™ *Tth* DNA Polymerase

Cat. Nos. TTH72100, TTH72250, TTH72500, TTH7201K, TTH7205K, and
TTH7225N

Connect with Epicentre on our blog (epicentral.blogspot.com),
Facebook (facebook.com/EpicentreBio), and Twitter ([@EpicentreBio](https://twitter.com/EpicentreBio)).

1. Introduction

MasterAmp™ *Tth* DNA Polymerase is a thermostable DNA Polymerase from *Thermus thermophilus*. The enzyme displays optimal activity at temperatures between 70°C and 74°C, and it is useful for performing the polymerase chain reaction (PCR)* or other reactions requiring DNA synthetic activity at high temperatures. *Tth* DNA polymerase has an intrinsic 5'→3' exonuclease activity, but lacks a 3'→5' (proofreading) nuclease activity. Furthermore, *Tth* has a potent Mn-dependent reverse transcriptase activity.¹ The enzyme is provided with MasterAmp *Tth* 20X PCR Buffer, MasterAmp 10X PCR Enhancer (with betaine)[†], and 25 mM MgCl₂, and MnSO₄ Solutions to allow optimization of individual template-primer pair combinations.

The presence of betaine (trimethyl glycine) in the MasterAmp 10X PCR Enhancer substantially improves the yield and specificity of amplification of many target sequences, especially those containing a high G+C content or secondary structure.²⁻⁶ Betaine, a natural osmoprotectant, lowers the melting temperature of G+C rich regions to a temperature more similar to A+T rich regions.⁷ This results in destabilization of double-stranded DNA which limits polymerase pausing,⁴ thereby increasing the yield of full-length product. In addition, betaine also may enhance PCR by protecting DNA polymerases from thermal denaturation.⁸ The effects of betaine seem to be independent of the polymerase used, though the concentration of betaine required for amplification varies with the target sequence.²⁻⁵

2. Product Specifications

Storage: Store only at -20°C in a freezer without a defrost cycle.

Storage Buffer: MasterAmp *Tth* DNA Polymerase is supplied in a 50% glycerol solution containing 50 mM Tris-HCl (pH 7.5), 100 mM NaCl, 0.1 mM EDTA, 0.5% Tween® 20, 0.5% NP-40, and 1 mM dithiothreitol.

Contaminating Activity Assays: MasterAmp *Tth* DNA Polymerase is free of detectable non-specific DNase and RNase activities.

Unit Definition: One unit converts 10 nmoles of deoxyribonucleoside triphosphates into acid-insoluble material in 30 minutes at 74°C using standard assay conditions.

Activity Assay: The activity assay is performed in a reaction containing 25 mM TAPS (pH 9.3), 50 mM KCl, 2.0 mM MgCl₂, 8.5 µg of activated calf thymus DNA, 0.2 mM of each dNTP, and 0.02-0.1 unit of enzyme.

MasterAmp *Tth* 20X PCR Buffer: 400 mM (NH₄)₂SO₄ and 1.0 M Tris-HCl (pH 9.0). Separate solutions of MasterAmp 10X PCR Enhancer, 25 mM MgCl₂, and 25 mM MnSO₄ are also provided to allow optimization of individual reactions.

Enzyme Characteristics:

5'→3' Exonuclease Activity Yes
 3'→5' Exonuclease Activity No
 Terminal Transferase Activity Yes
 Extension Rate (72°C) 2-4 kb/min
 Processivity 50-60 bases
 Half-life 10 min (97°C), 40 min (95°C)
 Reverse Transcriptase Activity (Mn²⁺ dependent) Yes

3. Kit Contents

Cat. #	Concentration	Quantity
The following include: MasterAmp <i>Tth</i> DNA Polymerase, MasterAmp <i>Tth</i> 20X PCR Buffer, MasterAmp 10X PCR Enhancer, 25 mM MgCl ₂ Solution, and MnSO ₄ Solution.		
TTH72100	@ 5 U/μl	100 U
TTH72250	@ 5 U/μl	250 U
TTH72500	@ 5 U/μl	500 U
TTH7201K	@ 5 U/μl	1,000 U
TTH7205K	@ 5 U/μl	5,000 U
MasterAmp <i>Tth</i> DNA Polymerase		
TTH7225N	@ 5 U/μl	250 U
MasterAmp 10X PCR Enhancer		
ME81201		1.5 ml
ME81205		5 ml
ME81210		10 ml
25 mM MgCl₂ Solution		
FB4250		25 ml

4. Related Products

The following products are also available:

- MasterAmp™ PCR Optimization Kits with Ammonium Sulfate
- MasterAmp™ PCR PreMixes
- MasterAmp™ AmpliTherm™, *Taq* and *Tfl* DNA Polymerases
- MasterAmp™ RT-PCR Kits
- MasterPure™ Nucleic Acid Purification Kits
- MasterAmp™ Buccal Swab DNA Extraction Kits
- BuccalAmp™ DNA Extraction Kits
- dNTP Solutions

5. General Considerations

1. **Template:** DNA or RNA prepared using standard isolation techniques are suitable substrates for amplification. Nevertheless, numerous compounds inhibit amplification including ionic detergents, some gel loading dyes, phenol, and hemin. When purifying templates from agarose gels, minimize exposure to UV irradiation to prevent formation of pyrimidine dimers. Assembly of reactions in a clean area or using positive displacement pipettors with aerosol-barrier tips will minimize the risk of contamination from extraneous DNA templates. The optimal amount of template for a single-copy gene is between 10⁴-10⁶ copies (e.g., approximately 0.1-10 ng of *E. coli* genomic DNA), though this may vary depending on the source and quality of the template.^{9,10}

- Primer Design:** Primers typically are 15-30 bases in length and contain approximately 50% G+C residues; the annealing temperatures of primer pairs should be nearly identical. Care must be taken to design primers that do not form hairpin loop structures or are self-complementary. The 5' end of a primer may contain bases that are not complementary with the template; however, the 3' end of the primer must be complementary with the template.
- Reaction Components:** DNA polymerases vary in their requirements for salt and pH, and therefore enzymatic activity will also vary depending on the reaction buffer. Mg^{2+} concentration is particularly critical for amplification of a specific target. Determine the optimal amount of Mg^{2+} for new template-primer pairs, or if the concentration of dNTPs and primers are altered. The ratio of primer to template is important for controlling the specificity and efficiency of amplification; an excess of primer ensures that the denatured template molecules bind to the primers instead of binding to the complementary DNA strand.¹¹ Nevertheless, too much primer may lead to the formation of nonspecific products or primer dimers.
- Choice of Enzyme:** Not all templates will be synthesized with equal efficiency by the various DNA polymerases. If amplification is not achieved with one enzyme, and after attempts to optimize fail, another enzyme should be chosen. Furthermore, use of a polymerase containing proofreading activity may be useful where fidelity or length of product is important.
- Cycling Parameters:** Many parameters influence both the specificity and efficiency of amplification including the temperature and duration of denaturation, annealing and elongation; ramp speed; and total cycle number. Amplification beyond $\sim 10^{12}$ molecules may also result in the appearance of nonspecific products; if the starting number of template molecules is 10^5 , 28-30 cycles will yield 10^{12} molecules.⁹ Additional variations such as use of a hot start¹² or touchdown/stepdown PCR^{13,14} can dramatically improve specificity and yield (see page 6 Troubleshooting Amplification Reactions).

6. Suggested PCR Protocol

Use the following protocol as a guideline for establishing the parameters necessary for amplification of experimental templates. The goals of the individual experiment (i.e., fidelity versus yield) will influence many aspects of the reaction, including the amount of various reaction components, primer design, cycling parameters and choice of DNA polymerase^{9,11,15} (see General Considerations). Assembly of amplification reactions in a clean area and using positive displacement pipettors with aerosol-barrier tips will minimize the risk of contamination from extraneous DNA templates.

A. Assemble the Reaction Mix

Preparation of a reaction mix simplifies the assembly of multiple amplification reactions. The volumes listed below are for one 50- μ l amplification reaction. Assemble an amount of mix corresponding to the total number of reactions. Extra mix may be required to offset losses caused by pipeting. We recommend duplicate reactions (+ and – MasterAmp 10X PCR Enhancer) for new template-primer pair combinations.

1. Thaw and thoroughly mix all of the reagents listed below before dispensing; place on ice.

Combine on ice, all of the following except for the DNA template (see Note below) in tubes labeled “+” and “-” MasterAmp Enhancer:

x	μl	sterile water	
2.5	μl	MasterAmp <i>Tth</i> 20X PCR Buffer	(1X final concentration)
2-10	μl	25 mM MgCl ₂	(1-5 mM final concentration)
0 or 15	μl	MasterAmp 10X PCR Enhancer	(3X final concentration)
4	μl	2.5 mM dNTP Mix	(200 μM final concentration each)
0.5-1.25	μl	20 μM Primer 1	(0.2-0.5 μM final concentration)
0.5-1.25	μl	20 μM Primer 2	(0.2-0.5 μM final concentration)
0.25	μl	MasterAmp <i>Tth</i> DNA Polymerase	(1.25 Units)
x	μl	DNA Template (see Note)	(1-1000 ng)
<hr/>			
50	μl	Total reaction volume	

Note: The DNA template is added directly to the individual PCR tubes and is not included in the tubes of reaction mix. (See step 2 below.)

B. Assemble the Reactions

1. Thoroughly mix the tube of reaction mix and aliquot into the individual PCR tubes.
2. Add the DNA template to the PCR tubes and mix.
3. Add mineral oil and centrifuge briefly if using a thermal cycler without a heated lid.
4. Program the thermal cycler following the recommendations provided by the manufacturer.

A suggested program is outlined below.

a) Initially denature the template at 92-95°C for 3-5 minutes.

b) Perform a 2- or 3-step cycling program, for 20-50 cycles as required:

Denature at 92°C-95°C for 0.5-1.0 minutes.

Anneal the primers at a temperature 5°C below the T_m of the primers for 0.5 minute.

Extend the annealed primers at 68°C-72°C for 1 minute for every kb of expected product.

c) A final extension may be performed at 68°C-72°C for 10-30 minutes to ensure full extension and to increase the likelihood of 3'-uncoded nucleotide addition (3'-terminal A overhang).

5. Place the tubes in the thermal cycler and begin cycling.
6. After amplification, the samples may be kept at 4°C overnight or frozen at -20°C.

7. Suggested RT-PCR Protocol

A. Assemble the Reaction Mix

Preparation of a reaction mix simplifies the assembly of multiple amplification reactions. The volumes listed below are for one 50- μ l amplification reaction. Assemble an amount of mix corresponding to the total number of reactions. Extra mix may be required to offset losses caused by pipeting. We recommend using 1X MasterAmp PCR Enhancer (final concentration) with new RT-PCR template-primer pair combinations. If successful amplification is not achieved, then perform subsequent reactions using either a final concentration of 0X or 2X MasterAmp PCR Enhancer. Note that the optimal concentrations of MasterAmp PCR Enhancer (0-2X), MnSO_4 (0.5-0.75 mM), MgCl_2 (1.5-3 mM), and dNTPs (200-400 μ M) may vary with each template.

1. Thaw and thoroughly mix all of the reagents listed below before dispensing; place on ice.

Combine on ice, all reagents in the order listed:

x μ l	sterile water	
2.5 μ l	MasterAmp <i>Tth</i> 20X PCR Buffer	(1X final concentration)
6 μ l	25 mM MgCl_2	(3 mM final concentration)
8 μ l	2.5 mM dNTP Mix	(400 μ M final concentration each)
5 μ l	MasterAmp 10X PCR Enhancer	(1X final concentration)
1 μ l	25 mM MnSO_4	(0.5 mM final concentration)
1.25 μ l	20 μ M Primer 1	(0.5 μ M final concentration)
1.25 μ l	20 μ M Primer 2	(0.5 μ M final concentration)
x μ l	RNA Template	(1-10 μ g)
0.5 μ l	MasterAmp <i>Tth</i> DNA Polymerase	(2.5 Units)
<hr/>		
50 μ l	Total reaction volume	

B. Assemble the Reactions

1. Thoroughly mix the tube of reaction mix and aliquot into the individual PCR tubes.
2. Add mineral oil and centrifuge briefly if using a thermal cycler without a heated lid.
3. For RNA templates up to 2 kb, synthesize the first strand by incubating the tubes at 60°C for 20 min. Add 10 minutes of incubation time for each additional kb.

If using a primer with a T_m of less than 60°C (e.g., oligo(dT)₁₈), incubate the reactions at an appropriate annealing temperature for 5 min before performing the synthesis at 60°C.

4. Program the thermal cycler following the recommendations provided by the manufacturer.

Perform a 2- or 3-step cycling program, for 30-50 cycles as required:

Denature at 92°C-95°C for 0.5-1.0 minutes.

Anneal the primers at a temperature 5°C below the T_m of the primers for 0.5 minute.

Extend the annealed primers at 68°C-72°C for 1 minute for every kb of expected product.

5. Place the tubes in the thermal cycler and begin cycling.
6. After amplification, the samples may be kept at 4°C overnight or frozen at -20°C.

8. Troubleshooting Amplification Reactions

Little or no amplification detected

- 1) **Lower annealing temperature.** Lower the annealing temperature in 2°C increments.
- 2) **Optimize Mg²⁺ concentration.** Perform reactions with varying concentrations of Mg²⁺ starting at 0.5 mM up to 5.0 mM, in 0.5 mM increments. Heat the vial of MgCl₂ at 70°C for 15 minutes and mix vigorously to resuspend any microcrystalline precipitates that may have formed.
- 3) **Perform hot start.**¹² The final assembly of amplification reactions at temperatures above the reaction annealing temperature improves both the yield and specificity. Assemble the reactions without Mg²⁺; subsequently place the reactions in a thermal cycler heated to >80°C, then add the appropriate amount of MgCl₂ and begin the cycling protocol. Alternatively, a modified hot start may be performed in which reactions assembled on ice are added directly to a thermal cycler pre-heated to 92-95°C.
- 4) **Perform Touchdown (TD)/Stepdown (SD) PCR.**^{9,13,14} TD or SD PCR consists of a series of cycles that are performed at decreasing annealing temperatures. This protocol favors amplification of the target at temperatures greater than or equal to the optimum annealing temperature, while enhancing target yield in later cycles at annealing temperatures below the T_m of the reaction. Perform the initial series of cycles at an annealing temperature a few degrees above the calculated T_m of the primers; annealing temperature is lowered by 1°C-4°C every other cycle to ~10°C below the calculated T_m. A hot start must be performed if using a TD or SD cycling protocol.
- 5) **Alter the concentration of MasterAmp PCR Enhancer in the reaction.** While 3X MasterAmp PCR Enhancer works well for many templates, concentrations of 1X, 2X, or 4X may work better.
- 6) **Increase initial template denaturation time or temperature.** Increase the temperature of initial denaturation up to 95°C. Increase the length of initial template denaturation up to 5 minutes. Alternatively, denature the template by heating at 72°C for 10 minutes in the presence of 50 mM NaOH before amplification.
- 7) **Increase number of cycles.** Perform additional cycles in increments of 5.
- 8) Vary reaction components. Vary the amount of DNA polymerase and primers. Try using an alternative DNA polymerase.
- 9) **Check template quantity and quality.** Check the concentration of template DNA by agarose gel electrophoresis or fluorimetry. Use 10⁴-10⁶ molecules of template for each reaction (e.g., up to ng amounts for cloned templates or µg amounts for genomic DNA).^{9,14} Organic extraction followed by ethanol precipitation may remove some inhibitors of amplification.
- 10) **Increase extension time.** Increase the extension time, generally 1 min for every kb of product.

Multiple products or a smear detected

- 1) **Decrease concentration of reaction components.** Check the concentration of template DNA by agarose gel electrophoresis or fluorimetry. Use 10^4 - 10^6 molecules of template for each reaction (e.g., up to ng amounts for cloned templates or μ g amounts for genomic DNA).^{9,14} Decrease the amount of magnesium, enzyme, and primer added to the reaction.
- 2) **Increase annealing temperature.** Increase the annealing temperature in 2°C increments.
- 3) **Perform hot start.**¹² The final assembly of amplification reactions at temperatures above the reaction annealing temperature improves both the yield and specificity. Assemble the reactions without Mg^{2+} ; subsequently place the reactions in a thermal cycler heated to $>80^\circ C$, then add the appropriate amount of $MgCl_2$ and begin the cycling protocol. Alternatively, a modified hot start may be performed in which reactions assembled on ice are added directly to a thermal cycler pre-heated to $92^\circ C$ - $95^\circ C$.
- 4) **Perform Touchdown (TD)/Stepdown (SD) PCR.**^{9,13,14} TD or SD PCR consists of a series of cycles that are performed at decreasing annealing temperatures. This protocol favors amplification of the target at temperatures greater than or equal to the optimum annealing temperature, while enhancing target yield in later cycles at annealing temperatures below the T_m of the reaction. Perform the initial series of cycles at an annealing temperature a few degrees above the calculated T_m of the primers; annealing temperature is lowered by $1^\circ C$ - $4^\circ C$ every other cycle to $\sim 10^\circ C$ below the calculated T_m .
A hot start must be performed if using a TD or SD cycling protocol.
- 5) **Alter the concentration of MasterAmp PCR Enhancer in the reaction.** While 3X MasterAmp PCR Enhancer works well for many templates, concentrations of 1X, 2X, or 4X may work better.
- 6) **Redesign primers.** Design primers that have higher annealing temperatures and do not form hairpin loops or primer dimers.
- 7) **Check primers for degradation.** Check by electrophoresis in a denaturing acrylamide gel.
- 8) **Decrease number of cycles.** Decrease number of cycles in increments of 5.

9. References

1. Myers, T.W. and Gelfand, D.H. (1991) *Biochem.* **30**, 7661.
2. Schanke, J.T. and Grunenwald, H.L. (1997) *Epicentre Forum* **4** (1), 2.
3. Grunenwald, H.L. and Schanke, J.T. (1997) *Epicentre Forum* **4** (1) 4.
4. Mytelka, D.S. and Chamberlin, M.J. (1996) *Nucl. Acids Res.* **24**, 2774.
5. Henke, W. *et al.*, (1997) *Nucl. Acids Res.* **25**, 3957.
6. Weissensteiner, T. and Lanchbury, J.S. (1996) *BioTechniques* **21**, 1102.
7. Rees, W.A. *et al.*, (1993) *Biochemistry* **32**, 137.
8. Santoro, M.M. *et al.*, (1992) *Biochemistry* **31**, 5278.
9. Roux, K.H. (1995) in: *PCR Primer*, Dieffenbach, C.W. and Dveksler, G.S. (eds.), CSH Press, New York, 53.
10. Kolmodin, L.A. and Williams, J.F. (1997) in: *PCR Cloning Protocols*, White, B.A. (ed.), Humana Press, Totowa, 6.

11. Cha, R.S. and Thilly, W.G. (1995) in: *PCR Primer*, Dieffenbach, C.W. and Dveksler, G.S. (eds.), CSH Press, New York, 37.
12. D'Aquila, R.T. et al., (1991) *Nucl. Acids Res.* **19**, 3749.
13. Don, R.H. et al., (1991) *Nucl. Acids Res.* **19**, 4008.
14. Hecker, K.H. and Roux, K.H. (1996) *BioTechniques* **20**, 478.
15. Kidd, K.K. and Ruano, G. (1995) in: *PCR 2: A Practical Approach*, McPherson, M.J. et al., (eds.), IRL Press, Oxford, 1.

AmpliTherm, BuccalAmp, MasterAmp, and MasterPure, and are trademarks of Epicentre, Madison, Wisconsin.

Tween is a registered trademark of ICI Americas, Wilmington, Delaware.

**Covered by German Patent No. DE4411588C1 and other pending patent applications in the U.S. and other countries assigned or exclusively licensed to Epicentre.*

**Notice to Purchasers: Limited License*

A license under US Patents 4,683,202, 4,683,195, and 4,965,188 or their foreign counterparts, owned by Hoffmann-La Roche Inc. and F. Hoffmann-La Roche Ltd. ("Roche"), has an up-front fee component and a running-royalty component. The purchase price of this product includes limited, nontransferable rights under the running-royalty component to use only this amount of the product to practice the Polymerase Chain Reaction ("PCR") and related processes described in said patents solely for the research and development activities of the purchaser when this product is used in conjunction with a thermal cycler whose use is covered by the up-front fee component. Rights to the up-front fee component must be obtained by the end user to have a complete license. These rights under the up-front fee component may be purchased from Perkin-Elmer or obtained by purchasing an Authorized Thermal Cycler. No right to perform or offer commercial services of any kind using PCR, including without limitation reporting the results of purchaser's activities or a fee or other commercial consideration, is hereby granted by implication or estoppel. Further information on purchasing licenses to practice the PCR Process may be obtained by contacting the Director of Licensing at The Perkin-Elmer Corporation, 850 Lincoln Centre Drive, Foster City, California 94404 or at Roche Molecular Systems, Inc., 1145 Atlantic Avenue, Alameda, California 94501 and outside the US at F. Hoffman-La Roche Ltd., Grenzacherstrasse 124, 4070 Basel, Switzerland.

Visit our technical blog: epicentral.blogspot.com