

## T4 DNA Ligase, Cloned

Cat. Nos. L0805H, L0810H, L0820H, LH805H, LH810H, and LH820H

Connect with Epicentre on our blog ([epicentral.blogspot.com](http://epicentral.blogspot.com)),  
Facebook ([facebook.com/EpicentreBio](https://facebook.com/EpicentreBio)), and Twitter ([@EpicentreBio](https://twitter.com/EpicentreBio)).

## 1. Introduction

T4 DNA Ligase is the most versatile and commonly used ligase for DNA cloning. This ATP-dependent enzyme covalently joins blunt or compatible cohesive ends, as well as nicks in double-stranded DNA. A 5'-phosphoryl group is required for ligation to a 3'-hydroxyl group. T4 DNA Ligase is available in 500-, 1,000-, and 2,000-Unit sizes at both standard (2 U/μl) and high (10 U/μl) concentrations. High-concentration T4 DNA Ligase is useful for obtaining maximum efficiency in blunt-end ligations. Both a 10X Reaction Buffer and 25-mM ATP Solution are provided with the enzyme.

For those interested in the shortest possible reaction times, Epicentre also offers Fast-Link™ DNA Ligation Kits. These kits are specially formulated to provide the fastest high-efficiency DNA ligations for routine and high-throughput DNA cloning. Cohesive-end ligations can be performed in as little as 5 minutes and blunt-end ligations in as little as 15 minutes with the Fast-Link Kits.

## 2. Product Specifications

**Storage:** Store only at -20°C in a freezer without a defrost cycle.

**Storage Buffer:** T4 DNA Ligase is supplied in a 50% glycerol solution containing 50 mM Tris-HCl (pH 7.5), 0.1 M NaCl, 0.1 mM EDTA, 1 mM dithiothreitol (DTT), and 0.1% Triton® X-100.

**Unit Definition:** One unit converts 1 nmol of pyrophosphate into Norit-adsorbable material in 20 minutes at 37°C.<sup>1</sup>

**Quality Control:** T4 DNA Ligase is function-tested in a reaction containing 33 mM Tris-acetate (pH 7.5), 66 mM potassium acetate, 10 mM magnesium acetate, 0.5 mM DTT, 1 mM ATP, 1 μg *Hind* III-cut lambda DNA, and varying amounts of enzyme.

**10X Reaction Buffer:** 330 mM Tris-acetate (pH 7.5), 660 mM potassium acetate, 100 mM magnesium acetate, and 5 mM DTT.

*ATP is not included in the 10X Reaction Buffer and must be added separately. For most applications, ATP should be added to the reaction to a final concentration of 0.5-1.0 mM.*

**Contaminating Activity Assays:** Cloned T4 DNA Ligase is free of detectable DNA exo- and endonuclease and RNase activities, as judged by gel electrophoresis following incubation of greater than 150 U of enzyme for 16 hours at 37°C with the following substrates: i) a 40-mer oligodeoxynucleotide with a 5'-phosphorylated end; ii) a 40-mer oligodeoxynucleotide with a 5'-hydroxyl end; iii) *Hind* III-digested bacteriophage lambda DNA; and iv) supercoiled pUC-19 DNA.

**Note on High-Concentration T4 DNA Ligase:** *At high concentrations, sufficient amounts of enzyme can bind to DNA, causing the DNA to migrate abnormally in an agarose gel and the DNA band to appear smeared. This phenomenon should not be mistaken as nuclease contamination, as the DNA is recovered intact if the Ligase is removed from the reaction. Excess T4 DNA Ligase can also decrease transformation efficiencies.<sup>2</sup> Excess enzyme may be removed by adding 0.1% SDS and Proteinase K (20 μg/ml) followed by incubation at 37°C for 30 minutes. The reaction is then heated to 70°C for 10 minutes, after which the DNA can be used in standard transformation protocols.*

### 3. Related Products

The following products are also available:

- Fast-Link™ DNA Ligation Kit
- Fast-Link™ DNA Ligation and Screening Kit
- Colony Fast-Screen™ Kits
- Ampligase™ Thermostable DNA Ligase
- T4 Polynucleotide Kinase
- T4 RNA Ligase
- T4 DNA Polymerase
- Mung Bean Nuclease
- End-It™ DNA End-Repair Kit

### 4. References

1. Weiss, B. *et al.*, (1968) *J. Biol. Chem.* **243**, 4543.
2. Michelsen, B.K. (1995) *Anal. Biochem.* **225**, 172.

*Triton* is a registered trademark of Rohm & Haas, Philadelphia, Pennsylvania.

*Ampligase* is a registered trademark, and *Fast-Link*, *Fast-Screen*, and *End-It* are trademarks of Epicentre, Madison, Wisconsin.

Visit our technical blog: [epicentral.blogspot.com](http://epicentral.blogspot.com)