

3. Parsons, L.M. *et al.* (1998) *Mol. Microbiol.* **28**, 571.
4. Goryshin, I.Y. *et al.* (2000) *Nature Biotechnol.* **18**, 97.
5. Berg, C.E. and Berg, D.E. (1996) in: *Escherichia coli and Salmonella: Cellular and Molecular Biology*, eds. Neidhardt, F.C. *et al.*, American Society for Microbiology, Washington, D.C., 2588.
6. Goryshin, I.Y. and Reznikoff, W.S. (1998) *J. Biol. Chem.* **273**, 7367.
7. Akerley, B.J. *et al.* (1998) *Proc. Natl. Acad. Sci. USA* **95**, 8927.
8. Griffin, T.J. *et al.* (1999) *Nucl. Acids Res.* **27**, 3859.
9. Gwinn, M.L. *et al.* (1997) *J. Bacteriol.* **179**, 7315.
10. Haapa, S. *et al.* (1999) *Nucl. Acids Res.* **27**, 2777.
11. Haniford, D.B. *et al.* (1991) *Cell* **64**, 171.
12. Chaconas, G. *et al.* (1996) *Curr. Biol.* **6**, 817.
13. Snapper, S.B. *et al.* (1990) *Mol. Microbiol.* **4**, 1911.
14. Pavelka, M.S. and Jacobs, W.R. (1996) *J. Bacteriol.* **178**, 6496.
15. Jacobs, W.R., Jr. *et al.* (1991) *Methods Enzymol.* **204**, 537.
16. Joyce, C.M. and Grindley, N.D.F. (1984) *J. Bacteriol.* **158**, 636.
17. Bardarov, S. *et al.* (1997) *Proc. Natl. Acad. Sci. USA* **94**, 10961.
18. Guilhot, C. *et al.* (1994) *J. Bacteriol.* **176**, 535.
19. Rubin, E.J. *et al.* (1999) *Proc. Natl. Acad. Sci. USA* **96**, 1645.
20. Machowski, E.E. *et al.* (2000), submitted to *Gene*.

EZ::TN™ <KAN-2>Tnp Transposome™

TSM99K2-F72 10 Reactions

Pre-formed Transposome containing the Tn903 kanamycin-resistance marker. The kit also includes 2 unlabeled primers.

EZ::TN™ <DHFR-1>Tnp Transposome™

TSM99D1-F72 10 Reactions \$ 350

Pre-formed Transposome containing dihydrofolate reductase gene (trimethoprim selection). The kit also includes 2 unlabeled primers.

Special Offer

Construct your own custom EZ::TN™ Transposon and make a Transposome™

Until 6/30/00 receive EZ::TN™ pMOD™<MCS> Transposon Construction Vector free when you purchase 10 Units of EZ::TN™ Transposase.

EZ::TN™ pMOD™<MCS> Transposon Construction Vector

MOD9201-F72 20 µg free with purchase of 10 Units of EZ::TN™ Transposase

EZ::TN™ Transposase

TNP92110-F72 10 Units

Visit our website at

www.epicentre.com/catalog/transposomics**Finish Your DNA Sequencing Project—****Complete Difficult Regions with SequiTherm™ EXCEL II DNA Sequencing Kits**

SequiTherm EXCEL II resolved a “2-year-old puzzle” due to “an unreadable hairpin loop” for the Washington University genome center.

- *Science*, Vol. 280, 8 May 1998, pg 816

Sequence your most difficult regions with SequiTherm™ EXCEL II DNA Sequencing Kits: hairpin loops, regions of high GC or AT content, areas of interstrand reannealing, and inverted or direct repeats. Using end labeled primers, SequiTherm™ EXCEL II DNA Sequencing Kits provide complete, clear sequencing data when all other methods fail.

Benefits

- Complete work on most any difficult gap to finish your sequencing project.
- High flexibility - you can perform both cycle and isothermal sequencing with one kit.
- Kits compatible with LI-COR® and ALF™ DNA Sequencers (both dye-primer and internal label) and with ABI™ automated sequencers (dye-primer).
- Reduced background signal.

G A T C



Figure. Sequencing through a GC-rich trinucleotide repeat. Supercoiled plasmid template containing (CGG)₂₃ was sequenced using the SequiTherm EXCEL II isothermal sequencing protocol.

SequiTherm EXCEL™ II DNA Sequencing Kit - LC For 25-41 cm gels on LI-COR® Sequencers

SE9101LC-F72 (100 Sequences)

SequiTherm EXCEL™ II DNA Sequencing Kit - LC For 66 cm gels on LI-COR® Sequencers

SE9202LC-F72 (100 Sequences)

SequiTherm EXCEL™ II Long-Read™ DNA Sequencing Kit - ALF™

For ALF DNA Sequencers

SE8301A-F72 (100 Sequences)

SequiTherm EXCEL™ II DNA Sequencing Kit (for manual cycle and non-cycle sequencing)

SEM79020-F72 20 Sequences

SEM79050-F72 50 Sequences

SEM79100-F72 100 Sequences